

PREPARATION OF HIGHLY PURIFIED PLANT DNA

phases are no longer evident. DNA will collect as white stringy mass. *Sample may be stored overnight at - 20°C at this point.*

9. Close and hook the end of a pasteur pipette and gently wind DNA onto end. Let liquid drain off and put into clean sterile tube containing 76% ethanol-10 mM ammonium acetate. Let stand for at least 20 minutes. DNA can be left on the pipette.

10. Remove DNA & pipette. Let liquid drain off (until almost dry & odor of ethanol is no longer evident) and

at atisopropylTj /F3 12 Tf 0 TcTD -0.0410il amla